Product number #10680

CII peptide ($CII_{259-273}$) – 100µg

- A Peptide for studies on T-Cell Responses Associated with Autoimmune Arthritis

Description

The CII peptide (CII259-273) contains the 259 to 273 sequence of rat CII. It is the non-galactosylated peptide that is ideal to serve as the control peptide for the Galactosylated CII peptide (CII259-273-GalHyl264; product numbers #10510 and #10550).

The Galactosylated CII peptide (CII₂₅₉₋₂₇₃-GalHyl₂₆₄; #10510 and #10550) is the dominant T-cell epitope in collagen induced arthritis (CIA) in mice¹. The Galactosylated CII peptide (CII₂₅₉₋₂₇₃-GalHyl₂₆₄) activates autoimmune T-cells when presented by the MHC II Aq protein². **The CII peptide** (#10680) **may be used as a control peptide** for the in vitro stimulus by the Galactosylated CII peptide (CII259-273-GalHyl264) in antigen-specific T-cell recall assays for quantitation of T cell activity 1,2,3,4,5 as determined by the selected readout by ELISA, ELISpot, flow cytometry or RNAseq, e.g.

Sequence

H2N-Gly-Ile-Ala-Gly-Phe-Lys-Gly-Glu-Gln-Gly-Pro-Lys-Gly-Glu-Thr-COOH

Peptide backbone sequence origin

Collagen alpha-1(II) chain, COL2A1, type II collagen, CII; UniProt: P05539.

Size

100 μg; á 20μl; (available also as 10 μg)

Supplied in

0.1M acetic acid

Molecular weight

1475.62 g/mol

Storage

It is recommended to store the peptide in aliquots at -20°C, or preferably at -80°C, unless stability data would suggest otherwise. If the peptide is stored less than 48 hours before use, the peptide may be stored at 5°C (or room temperature). Avoid repeated thawing and freezing procedures.

Recommended dilution

It is recommended that the user determines the optimal dilution for their application. For in vitro T cell stimulation, a concentration of 10 $\mu g/ml^5$ and a concentration of 20 $\mu g/ml$ for IL-17 ELISpot assays⁴ has been used.

For Research Use Only. Not for use in diagnostic procedures. Not for resale without express authorization.

References

- Michaëlsson E, Andersson M, Engström A, Holmdahl R. Identification of an immunodominant type-II collagen peptide recognized by T cells in H-2q mice: self tolerance at the level of determinant selection. Eur J Immunol. 1992 Jul;22(7):1819-25. PMID: <u>1378019</u>
- 2. Andersson IE, Andersson CD, Batsalova T, Dzhambazov B, Holmdahl R, Kihlberg J, Linusson A. Design of glycopeptides used to investigate class II MHC binding and T-cell responses associated with autoimmune arthritis. PLoS One. 2011 Mar 15;6(3):e17881. PMID: 21423632.
- 3. Lindgren C, Andersson IE, Berg L, Dobritzsch D, Ge C, Haag S, Uciechowska U, Holmdahl R, Kihlberg J, Linusson A. Hydroxyethylene isosteres introduced in type II collagen fragments substantially alter the structure and dynamics of class II MHC A(q)/glycopeptide complexes. Org Biomol Chem. 2015 Jun 14;13(22):6203-16.. PMID: 25960177.
- 4. Klocke K, Sakaguchi S, Holmdahl R, Wing K. Induction of autoimmune disease by deletion of CTLA-4 in mice in adulthood. Proc Natl Acad Sci U S A. 2016 Apr 26;113(17):E2383-92. PMID: 27071130.
- Merky P, Batsalova T, Bockermann R, Dzhambazov B, Sehnert B, Burkhardt H, Bäcklund J. Visualization and phenotyping of proinflammatory antigen-specific T cells during collagen-induced arthritis in a mouse with a fixed collagen type II-specific transgenic Tcell receptor β-chain. Arthritis Res Ther. 2010;12(4):R155. PMID: 20682070

Vacara AB

Nanna Svartz väg 4, 171 65, Solna, Sweden

Tel: +46 70 765 6638 Email: info@vacara.se

